

Enteric coatings with EUDRAGIT[®] L/S from organic solutions. Tablets coated by fluid-bed process

Spraying process for the manufacture a coating of colourless enteric sealing coats with EUDRAGIT[®] L 12.5 on tablets.

Operating method

The coating suspension is sprayed onto the fluidized tablets, which are prewarmed to about 30 °C, by means of built-in air nozzles (bottom-spray method). Spray rate, inlet air quantity and inlet air temperature are adjusted in such a way that spraying can be performed continuously. During the process, the tablets should be maintained at a temperature of 25 to 30 °C and be able to flow freely.

Twinning, i.e. sticking together of tablets, can be avoided by adding suitable glidants (talc, magnesium stearate, kaolin, Syloid[®]) to the EUDRAGIT[®] L 12.5 spray suspension.

If twinning does occur, spraying must be interrupted until the tablets are dry and once more able to flow freely. Subsequently, processing may be continued at a reduced spray rate.

The following formulation gives the polymer and excipient quantities required for coating 1 kg of medium-sized tablets (diameter 11 mm, weight 640 mg) at a polymer weight of 8.1 mg/cm².

The solids are combined with the acetone and EUDRAGIT[®] L 12.5 and are then finely dispersed in a doublecone mill.

EUDRAGIT[®] L 12.5 is also easy to prepare from the solvent free solid EUDRAGIT[®] L 100. To this end, 13 parts by weight of EUDRAGIT[®] L 100 (= 12.5% dry polymer substance) are dissolved in a mixture of 82 wt. parts isopropyl alcohol and 5 wt. parts water at room temperature with stirring.

Typical formulation from organic solution

	Parts by weight
EUDRAGIT[®] L 12.5	440 g
Plasticizer	5 g
Talc	15 g
Elcema [®] P 050	5 g
Acetone	535 g
	à 1,000 g
Solids content:	8.0%
Content in dry polymer substance:	5.5%

This gives a polymer (55 g from 440 g EUDRAGIT[®] L 12.5) weight of $\Gamma = 8.1 \text{ mg/cm}^2$ by a tablet surface of $S = 432 \text{ mm}^2$.

Where delayed release in the intestine is required in addition to resistance to gastric fluid, EUDRAGIT[®] L can be partially or completely replaced with EUDRAGIT[®] S.

Dosage forms with pH-independent delayed release of the active ingredient are obtained by using EUDRAGIT[®] RL/RS as the film formers. If organic solvents must be avoided altogether, the aqueous dispersion EUDRAGIT[®] L 30 D-55 is a suitable alternative.

Operating data

Example

Technical data

Coating unit

EUDRAGIT® L/S

Aeromatic STREA-1 single-chamber fluidized-bed coater, bottom-spray

Feed pump

peristaltic pump with silicone tube, internal dia. 3 mm

Spray system

air spray gun, nozzle dia. 1.1 mm (round spray)

distance nozzle/tablets

120 mm

Coating data

Tablets

active cores, 11 mm in dia., 7 mm in height, 640 g in weight; hardness 70 N, disintegration time in demineralized H₂O: 30 sec

Batch size

1 kg

Spray suspension

620 g, corresponding to 50 g solids or 5 wt.-% on tablet quantity, 34 g polymer equivalent to 3.4 wt.-% on tablet quantity, ratio 2:1 to other excipients

Process data

Duration

Preheating

2 min

Spraying

21 min

Drying

3 min

Inlet air quantity

-

2 m³/min

1.6 m³/min

Inlet air temperature

37 °C

36 °C

36 °C

Air inlet setting

5

13

11

Outlet air temperature

30 °C

29 °C

30 °C

Tablet temperature

30 °C

29 °C

30 °C

Pump speed

-

10 rpm

-

Atomizing air pressure

2 bar

2 bar

2 bar

Spray rate

-

30 g/min

-

Other process data

Spray rate

30 g suspension/min/kg product = 2.38 g solids/min/kg product

Evaporation rate

27.1 g/min/kg product

Polymer quantity

5.0 mg/cm²

Spraying process

continuous

Total process time

26 min

Post-drying

2 hrs on trays, in drying cabinet at 40 °C

Results

Appearance

uniform coating of subdued gloss

Gastroresistance

for 3 hours (BP), disintegration in intestinal fluid (BP): max. 45 min

Our technical advice on the applications of our products is given without obligation. The buyer is responsible for the use and processing of our products and is also liable for observing any third-party rights. Technical data concerning our products are typical values. Subject to alteration.

® = registered trademark

EUDRAGIT = reg. Trademark of Röhm GmbH & Co. KG, Darmstadt, Germany